ANGELINA COLLEGE SAFETY-SHOWER/EYE WASH INSPECTION INSTRUCTIONS

NO. El-03

TITLE: Safety Shower/ Eye Wash Inspection Instructions

6.3.a

PAGE 1 of 1 **ISSUED**: 10/15

REVISED: 6/27

NOTES:

- Eye washes must be tested and inspected weekly.
- An Eye Wash Inspection Tag will be attached to each eye wash.
- Outlet heads (lids covering where water flows from) should be kept closed when not in use

INSTRUCTIONS:

- Run the eye wash for 2-3 minutes, ensuring:
 - o there is sufficient water flow
 - o the hands-free mechanism is functioning
- Initial the appropriate box on the attached Inspection Tag to document a passing inspection
- If inspection fails, notify all users and call Campus Services: 633-5280 for repairs
 - Should an exposure occur, flush the affected eye(s) for 15 minutes.
 - To ensure adequate flushing, hold eyelid(s) open and roll the eyeball.

INSPECTION TAGS			
Date of Inspection	Initials of Person Testing	Date of Inspection	Initials of Person Testing